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## Instruction for Use TPP Tissue Culture Test Plate

*TPP tissue culture plates are designed for the cultivation and growth of cells as well as for photometric measurements and cell-based assays. The growth areas - but not the sidewall of the wells - are opto-mechanically activated for optimal cell adhesion and growth.*

*The tissue culture test plate is intended for single use only.*

### Note:

*Follow national regulations when handling biological material, use appropriate protective clothing*

*Take note of the regulations of aseptic working during the work process.*

### Instructions:

- *Check the packaging material and the product for intactness. Use only faultless material.*
- *Check the expiry date (EXP) on label and packaging. Only use products with valid EXP.*
- *Open the packaging in a sterile environment and remove a complete system (lid with plate) for use.*
- *Open the plate and fill it with the medium and inoculum according to the usual laboratory routine.*
- *Cover the wells with the supplied lid to reduce the risk of cross-contamination.*
- *Cultivate and harvest the cells according to the usual laboratory routine.*
- *In order to achieve uniform growth on the surface, the following aspects must be considered:*
  - *Resuspend the cells to avoid cell clumps. (Uneven growth)*
  - *Avoid foaming during resuspension creating holes in the cell lawn.*
  - *Low medium volume leads to the formation of minuscule cells, which leads to an uneven growth of the cells. (many cells on the edge, few in the middle).*

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#### **Advice and Recommendations:**

- *The stacking rim enables safe stacking of several plates, also in combination with other TPP plates of different well number.*
- *For a homogenous cultivation of the cells, levelling of the incubator and continuous supply of medium is essential.*
- *Avoid and/or eliminate vibration sources near the incubator to ensure uniform cell growth.*

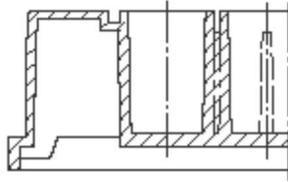
**ELISA:** The TPP plates feature are **not** suited for the use in ELISA tests.

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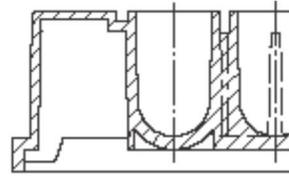
**Additional:**

Instructions for use, chemical resistance lists, and quality certificates of the respective products are available for download on the homepage [www.tpp.ch](http://www.tpp.ch).